

Instructions for Dino-Lite AM-312/412 for Television

Connecting to a television:

It is recommended using an LCD television or Plasma television for better image results. Older generation televisions with video input features will work also. For televisions without direct video input jack, an external device such as a VCR or DVD player with video input can be utilized.

Follow this procedure to connect model AM-312/412 to a television.

IMPORTANT: *The television must have an available video input jack to complete this procedure.*

1. Connect the composite video cable (yellow) to an available **VIDEO** input jack (yellow) on the television (Figure1).
2. Connect the female power cable end to the supplied male AC power adapter (Figure 2).
3. Plug in the AC power adapter into a power outlet. (Figure 3)
4. Turn on the television.
5. Tune the television to the appropriate video input channel to receive the signal from the digital microscope.

NOTE: *If necessary, consult the instruction manual of your television for information about tuning the TV to the correct video input channel.*

How to use:

Simply turn the adjustment dial left or right to adjust focus and magnification.

Included in package

- Dino-Lite digital microscope AM312/412
- AC power adapter (110V)
- Cradle stand



Options:

- Light Diffuser (sold separately) for highly reflective surfaces

Figure 1



Figure 2



Figure 3

